HAMILT®N

Intelligent Sensors

Arc Gives You the Power to Connect



Arc: The Power to Connect **RETHINK SENSOR MANAGEMENT BY ELIMINATING THE TRANSMITTER**

Arc seamlessly connects with existing R&D, PD, and production infrastructure and reduces startup costs and integration footprint on new projects and expansions. ArcAir software reduces operational costs with:











LABORATORY CALIBRATION

PREDICTIVE DIAGNOSTICS



USER ROLES AND PROCESS ASSIGNMENT



ANATOMY OF AN ARC SENSOR

Arc's integrated microtransmitter amplifies sensor signal for direct connection to the control system. Sensor configuration and calibration occur via USB or wireless Bluetooth®



Arc sensors save space and cost with their integrated microtransmitter. Calibration and diagnostic data is saved within the sensor, allowing for more reliable in-lab calibration.

Hardwired Process Control Data

Arc sensors provide a robust connection directly to the Process Control System. A wide array of analog and digital communication protocols allows effortless integration into existing systems:

• 4-20 mA • Profibus Modbus Foundation Field Bus

Arc Wi Adapter (optional)

Arc Wi Adapters enable wireless communication from the sensors to the ArcAir application without interrupting the hardwired signal. An LED in the adapter provides visual indication of sensor status.

Status Lights

 Active BT Connection
 Sensor Warning Normal Operation Sensor Errors



Wireless Configuration, **Calibration, and Diagnostics**

Arc Wi provides wireless communication to any Bluetooth 4.0 enabled device. Configure and calibrate sensors from the convenience of a mobile phone, tablet, or computer. Sensor diagnostics are also transmitted wirelessly, providing the ability to troubleshoot sensors from the palm of your hand.



| 2:05 PM | ······································ |
|------------------------|--|
| € \9 | Search |
| VCD | Incyte Arc |
| 5.86e6 cells/ml | 23.74°C |
| • pH | 238999-5241-1002 |
| 5.39pH | 23.09°C |
| • pH | 242120_0002627 |
| 7.27pH | 23.43°C |
| • DO | 242163-5187 LW |
| 97.95%-sat | 26.92C |
| DO | 243666-111-1178 |
| 92.76%-sat | 26.90C |
| • pH | PolilytePlus TS2 |
| 2.72pH | 22.79C |
| • ORP | 243061-erx |
| 93.20mV | 296.69K |
| DO | 243110-1078 |
| 0.00%-vol | 602.22C |
| Cond 0.01uS/cm | 243590-111-1258 22.98°C |
| • VCD | Incyte Arc |
| 0.00 | 23.58°C |
| Cond | RO Water #1 |
| 0.11uS/cm | 24.04°C |
| Cond | 242159-3357 |
| 0.01uS/cm | 22.68°C |
| Cond | 242159-3946 LW 🗙 |
| ? ⊵ | |

ArcAir for PC and Mobile ONE TOOL FOR SENSOR MANAGEMENT AND DOCUMENTATION

ArcAir is intuitive software designed to work the way you do. Monitor all sensors in your lab from an easy-to-use dashboard with one-click access to sensor configuration, calibration, and verification. Maintenance actions are automatically recorded, and reports are generated from a centralized database.

ArcAir Software Packages


| Functionalities | ArcAir Basic | ArcAir Advanced |
|--------------------------------------|--|--|
| Sensor Status | Image: A second s | Image: A second s |
| View and Sort Dashboard | Image: A second s | Image: A set of the set of the |
| View Experiments | Image: A second s | × |
| Create Experiments | Image: A second s | Image: A second s |
| Multi-language | Image: A set of the set of the | × |
| Configure / Calibrate / Verify | Image: A set of the set of the | Image: A second s |
| Firmware Updates | Image: A second s | × |
| Audit Trail with Comment Function | | Image: A set of the set of the |
| GMP Documentation Support | | × |
| User's Roles and Password Management | | × |
| Encrypted Database Backup and Change | | × |
| Reports Electronic Signatures | | × |

Mobile app is available through application stores such as App Store® or Google Play[™], with in-app upgrades for Basic and Advanced. Lite version is mobile app only. Basic and Advanced are available for PC. PC software available at hamiltoncompany.com/ArcAirSoftware.



Connection Flexibility

ArcAir is compatible with Android[™], iOS[®], and Windows[®]. Connect via USB or Bluetooth 4.0.

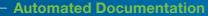

Eliminate manual reporting for more detailed, error-free documentation in less time.

- Print documentation for sign-off or save as PDF
- GMP-compatible reporting
- Mobile data synced to centralized database
- Searchable report history



-

-

-

atting a

12.00

Quality reduces INT & Quality Sector Biol 71 & Manufacture 2010 1



Configuration and Calibration

Easy-to-follow wizards and quick views simplify everyday tasks.

- Mobile barcode scanning for buffer lot tracking
- Graphical calibration, verification, and communication validation wizards
- Quickly view sensor health and error messages on one screen
- Cloning of sensor configurations

Common Interface for Mobile and PC

Arc sensor dashboard is a concise view of all sensors in the lab. Create sensor views by grouping and filtering. Access any sensor with one click from the dashboard.

Redefine Your Sensor Workflow

HASSLE-FREE SENSOR HANDLING

Arc breaks the reliance on at-line calibrations during start-up, takes the guesswork out of qualifying a used sensor for the next run, and eliminates the burden of manual documentation.

Grab and Go

Precalibrated sensors are waiting on the shelf for installation and sterilization with no additional at-line calibration.

Batch Calibration

Arc microtransmitters store calibration data in the sensor so you can calibrate on your time.

A dedicated batch calibration station increases efficiency and accuracy of calibrations.



ArcAir Dashboard

Track the status of all sensors in the lab on a single screen. Colored indicators enable quick identification of sensors needing attention.

Automatic generation of calibration and validation reports removes the risk of

transcription errors. Electronic signatures

enable a completely paperless workflow.

Historic sensor data can be used for troubleshooting or for predicting the useful life of a sensor in a specific process. Part 11

Dead Sensors

Sensor health indicators, operating hours, and sterilization counters indicate if the sensor should be discarded or if it is adequate for another run.

Arc Database

ArcAir stores all calibration, validation, and configuration information in an encrypted database. All user entries and actions are logged from all devices in an automatically generated audit trail.

Compliance Ready

Ready for compliance with FDA CFR 21 Part 11 and EudraLex Volume 4 Annex 11.



Mobile Friendly

Wirelessly receive diagnostic information from the sensors without interrupting the process. Quickly sync information to the database for reporting and data analysis.

Arc for GMP Facilities **INCREASE PROCESS SECURITY** AND AUTOMATE DOCUMENTATION

Traditional transmitters do not support a modern security model, leaving measuring points susceptible to uncontrolled user access. ArcAir Advanced accommodates the security and documentation needs of GMP facilities with a centralized database for distribution of user roles and storage of sensor history.



User Roles and Process Assignment

ArcAir software enables administrators to create optional user accounts and access levels for enhanced security. After login, users can only access assigned sensor groups and tasks within their role.

| Role: | Role Rights: | CFR 21 Part 11 and EU Annex 11 Compliance: | | Sensor Process Group Assignment Example: | | | Arc Database | |
|-------------|---|---|--|---|------------|------------|--|--------------------|
| | | ArcAir Software Settings | Electronic Signature For Report Creation/Approval | TANK 1 | TANK 2 | TANK 3 | | |
| ADMIN | AdministratorSensor SettingsProcess SettingsFirmware UpdatesCalibrationUser ManagementAudit Trail ViewCommunication Validation | | | ~ | ~ | ~ | Audit Trail | |
| P TECH 1 | Calibration Technician • Communication Validation • Process Setting • Audit Trail View | \bigcirc | \checkmark | \bigcirc | | \bigcirc | All user's GMP relevant activitie performed with mobile and desl are automatical | es Arc sktop |
| C TECH 2 | Production Technician Communication Validation Audit Trail View Process Settings | \bigcirc | | | \bigcirc | | logged in the au | |
| QM | Quality Manager Audit Trail View | \bigcirc | | \bigcirc | \bigcirc | \bigcirc | | |

الرق الرق

CONFIGURATION REPORT

> CALIBRATION REPORT

VERIFICATION REPORT

COMMUNICATION

REPORT

Process Documentation

ArcAir automatically generates and archives pre- and post- process documentation.

User and Sensor Configuration

Configure new sensors with ArcAir

Sensor Calibration

Calibrate sensors under controlled laboratory conditions

Communication Validation

Verify signal accuracy to the PCS

Process

Capture and export process measurement data

Sensor Verification

Verify sensor performance after process

Small Footprint, Big Impact DEMAND HIGHER VALUE FROM YOUR SYSTEM DESIGN

Modern equipment manufacturers implement Arc sensors to save space, improve usability, and expedite validation, ultimately reducing project installation and operating costs.

Break from the Bull

Shrink your panel and minimize failure points by eliminating bulky transmitters. Wiring is simplified to reduce assembly and troubleshooting costs.

Powered by Arc -

Accelerate feature-rich controls integration with direct access via Modbus RTU protocol. Leverage built-in algorithms for simple implementation of complex tasks like sensor health indication, error and warning descriptions, and calibration and validation variables. TCD

VCD

More Robust IQ/OQ/PQ

Skid builders capture performance data using the ArcAir Experiments tool for enriched Factory Acceptance Testing (FAT) and Functional Performance Testing (FPT). Arc's automated reports fast-track Turn Over Package (TOP) creation for rapid on-site validation.

Green Means Go

Quickly identify sensor status from a distance with colored LEDs. Use wireless diagnostics and calibration through a mobile or PC to remotely troubleshoot when needed.



Ò

0

0

0

©2022 Hamilton Company. All rights reserved.

Bluetooth is a registered trademark of Bluetooth SIG, Inc. App Store, iOS, and their logos are registered trademarks of Apple Inc. in the United States and other countries. Android, Google Play, and their logos are trademarks of Google, Inc. in the US and/or other countries. Windows and their logos are registered trademarks of Microsoft Corporation in the United States and other countries. All other trademarks are owned and/or registered by Hamilton Company in the U.S. and/or other countries. REF 691069/07 - - 08/2022

Ò



Web: www.hamiltoncompany.com USA: 800-648-5950 Europe: +41-58-610-10-10

Hamilton Americas & Pacific Rim

Hamilton Company Inc. 4970 Energy Way Reno, Nevada 8950<u>2 USA</u> Tel: +1-775-858-3000 Fax: +1-775-856-7259 sales@hamiltoncompany.com Hamilton Europe, Asia & Africa Hamilton Bonaduz A.G. Via Crusch 8 CH-7402 Bonaduz, Switzerland Tel: +41-58-610-10-10 contact.pa.ch@hamilton.ch

To find a representative in your area, please visit hamiltoncompany.com/contacts.