



SoftGrip™ Multi-Channel Pipettes

User Manual



Table of Contents

Welcome	1
Pipette Warranty	1
Hamilton SoftGrip™ Pipettes	3
Volume Setting	4
Recommended Volume Ranges	4
Pipette Operation	5
Adjusting for Left-Handed Operation	5
Pre-Rinsing Tips	5
Pipette Calibration	5
To Calibrate	6
Calibration Label and Key	7
Performance Specifications	7
Optimizing Performance	8
Troubleshooting	8

Cleaning	9
Autoclaving	9
Pipette Parts Diagrams	9
Replacement Parts Table	9
Replacement Parts	10
Pipette Recertification and Repairs	11
Return of Goods	12
Hamilton Precision Pipette Tips	12

Welcome

Thank you for choosing a SoftGrip™ pipette from the Hamilton Company.

Before using your new pipette, please review the important information in this manual on pipette features, basic operation, calibration, optimizing performance, maintenance, and troubleshooting.

For French, German, and Spanish translations of this manual, please visit www.hamiltoncompany.com/pipettes. For sales and support in countries outside the U.S., please contact your local dealer.

For additional information in the U.S., call Hamilton Sales and Support at 1-888-525-2123 and visit our website at www.hamiltoncompany.com/pipettes.

Pipette Warranty

Hamilton Company unconditionally guarantees its products to be free of defects in materials and workmanship for one year from date of delivery. Any product that fails due to such defects will be repaired or replaced at our discretion without cost, provided the device is returned with an explanation. It is the responsibility of the purchaser to determine the suitability of application and material compatibility of the products based on the published specifications of the products. The specified accuracy and precision performance of Hamilton pipettes is guaranteed only when used with Hamilton Precision Pipette Tips.

SoftGrip pipettes are protected by U.S. Patent Number 5,983,733 with other patents pending.

User Manual

SoftGrip Multi-Channel Pipettes

Hamilton SoftGrip™ Pipettes

Volume Setting

Recommended Volume Ranges

Pipette Operation

Adjusting for Left-Handed Operation

Pre-Rinsing Tips

Pipette Calibration

To Calibrate

Calibration Label and Key

Performance Specifications

Optimizing Performance

Troubleshooting

Cleaning

Autoclaving

Pipette Parts Diagrams

Replacement Parts Table

Replacement Parts

Pipette Recertification and Repairs

Return of Goods

Hamilton Precision Pipette Tips


Hamilton SoftGrip™ Pipettes

Hamilton Company is the leader in precision fluid measurement products. We extended our world-renowned syringe expertise to the SoftGrip Manual Pipettes specifically designed to deliver maximum accuracy and minimize hand injuries. The results set the standard for precision, comfort, and value.

SoftGrip multi-channel pipettes are continuously adjustable volume pipettes and available as 8- or 12-channel models for pipetting into 96-well plates. The pipettes feature a soft, nonslip handle for a relaxed grip; lighter plunger action for more comfort; and a separate tip ejector to reduce hand motion. The curved hilt supports the pipette for finer control while pipetting. The easily activated tip ejector lever releases all tips with a single stroke. The ejector lever rotates to a comfortable setting for left- or right-handed use with a simple turn of the lower body adjustment ring and tip ejector.

Each channel has a separate piston and seal. The hardened, stainless steel pistons are precision ground and polished to ensure accurate performance, and provide exceptional piston and seal life.

SoftGrip pipettes operate on the air displacement principle and use disposable tips. Every pipette is calibrated at the factory with standards traceable to N.I.S.T. and shipped with a Certificate of Calibration documenting the pipette's accuracy and precision.

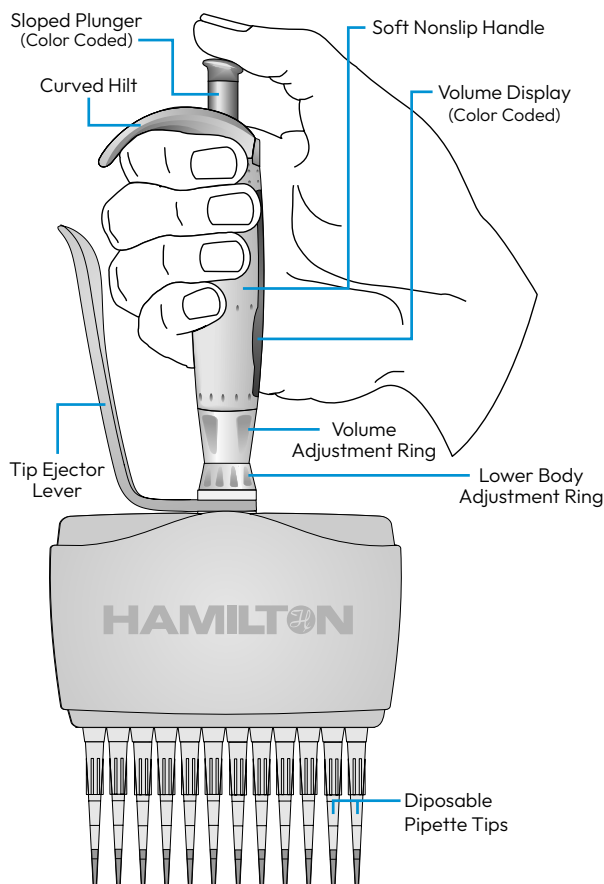

Volume Setting

The isolated volume adjustment ring prevents accidental volume changes. To set the volume, hold the pipette vertical with the volume display in front of you and turn the volume adjustment ring to the desired volume. Set the desired volume to within the pipette's volume range as given below.

For the greatest attainable accuracy, always turn the adjustment ring down to the desired volume. If you are starting at a setting lower than the one desired, turn the adjustment ring past that volume and then turn back to the desired volume.

When setting volumes, start your settings from the top digit on the volume setting display. Full units are expressed in black digits. Incremental units, the values to right of the decimal point, are expressed in red digits and appear below the full units. The graduated scale identifies the least significant digit.



Note: For the 50 μL pipettes, the red digits on the scale reading are to the right of the decimal point.

For the 300 μL pipettes, only the last digit, obtained from the graduated scale, is to the right of the decimal point.

For the best accuracy and precision, use your pipette only within the recommended volume range of 10-100% of the specific pipette's total volume range.

Recommended Volume Ranges

Pipette Volume	50 μL	300 μL
Pipette Color	Sandstone	Brick Red
Recommended Volume Range	5-50 μL	30-300 μL
100% Volume Setting	50.00 μL	300.0 μL
Reading at 100% Volume		
10% Volume Setting	5.00 μL	30.0 μL
Reading at 10% Setting		
Smallest Increment of the Scale	0.02 μL	0.2 μL

Pipette Operation

1. Press disposable tips firmly onto the pipette's lower body assembly. To assure an airtight seal, gently rock the handle from side to side before lifting the tips from the rack.
2. Depress the plunger button to the first stop. Hold the pipette vertically and immerse the tips 2 to 4 mm into the sample below the meniscus.
3. Slowly release the plunger button to aspirate sample.
4. Withdraw the tips from the sample, touching-off on the side of the container.
5. To dispense, place the tips against the inside wall of the receiving vessel and depress the plunger button to the first stop. Pause and then depress the plunger button to the second or blowout stop. Hold the plunger button down, sliding the tips out along the inside of the receiving vessel.
6. Release the plunger button. With the disposable tips directed into a safe receptacle, depress the tip ejector to dispose of the tips.

Adjusting for Left-Handed Operation

The pipette is delivered ready for right-handed operation. Both the tip ejector and lower body can be easily rotated for left-handed operation.

1. The tip ejector lever can be rotated a full 360° relative to the handle by gently pushing the lever away from the handle. While the handle assembly is lifted, rotate the lever to the desired position and then release the lever.
2. The entire lower body assembly can be rotated a full 360° relative to the handle by loosening the lower body adjustment ring. Rotate the lower body to the desired position and retighten the lower body adjustment ring.

Pre-Rinsing Tips

Pre-rinsing pipette tips is recommended for all pipetting operations. To pre-rinse, aspirate and dispense the sample one to three cycles. On the last rinse, press the plunger past the blowout stop to clear the tip of any sample. Make sure there are no micro-droplets in the tip before aspirating the first sample. If micro-droplets remain in the tip, replace the tip and pre-rinse again.

Pipette Calibration

Hamilton SoftGrip pipettes can be gravimetrically calibrated by the user. Each pipette comes with its own calibration key so that the pipette can be adjusted to accommodate specific fluid viscosities and environmental factors such as temperature, humidity, altitude, etc.

Use a 5- or 6-place balance depending on the volume being measured. A 5-place balance should be accurate to within ± 0.01 mg and a 6-place balance should be accurate to within a ± 0.001 mg. All items used to perform the calibration must be allowed to equilibrate to room temperature (20–25°C, constant to $\pm 0.5^\circ\text{C}$).


Each time you recalibrate the pipette, initial and date a new calibration label and replace the current label. Also, in accordance with your organization's instrument operation practices and standards, we recommend following good laboratory practices by calibrating your pipette on a regular basis, i.e., three month intervals, using the following procedure, or a known standard method, such as DIN 12650. Depending on the frequency of use and the liquids measured, other time intervals may be more appropriate.

To Calibrate

1. Attach a disposable tip to a chosen channel (a middle channel is recommend) and pre-rinse the tip.
2. Set the desired or the lowest volume at which the pipette will be used.
3. Make 10 measurements of distilled water per tip, recording the weights. Adjust for environmental factors, e.g., temperature, humidity, and altitude.
4. Calculate the mean and standard deviation.
5. If necessary, adjust the volume. However, adjust the volume only after confirming the error is not due to other causes. Refer to the sections on Optimizing Performance and Troubleshooting. Adjusting the volume will affect all channels.
 - a. Remove the calibration label and locate the calibration reference mark.
 - b. Remove the plunger cap by pulling the plunger button off the pipette.
 - c. Inset the calibration key, turning the calibration key clockwise to increase volume and counter-clockwise to decrease volume. Each vertical line on the calibration key is approximately 0.2% of the pipette volume.
 - d. Remove the calibration key and replace the plunger cap.
Repeat the gravimetric procedure until results are satisfactory.
6. Initial and date a new calibration label, placing the label on the pipette hilt.

The pipette can be recalibrated for liquids with a density other than water. Follow the calibration procedure above with the pipette set at the required volume. Use the liquid to be sampled. With this type of calibration adjustment, the pipette is calibrated only for the liquid used and the volume tested. The pipette should be labeled as such, and should not be used for other sampling applications unless it is recalibrated.

Calibration Label and Key

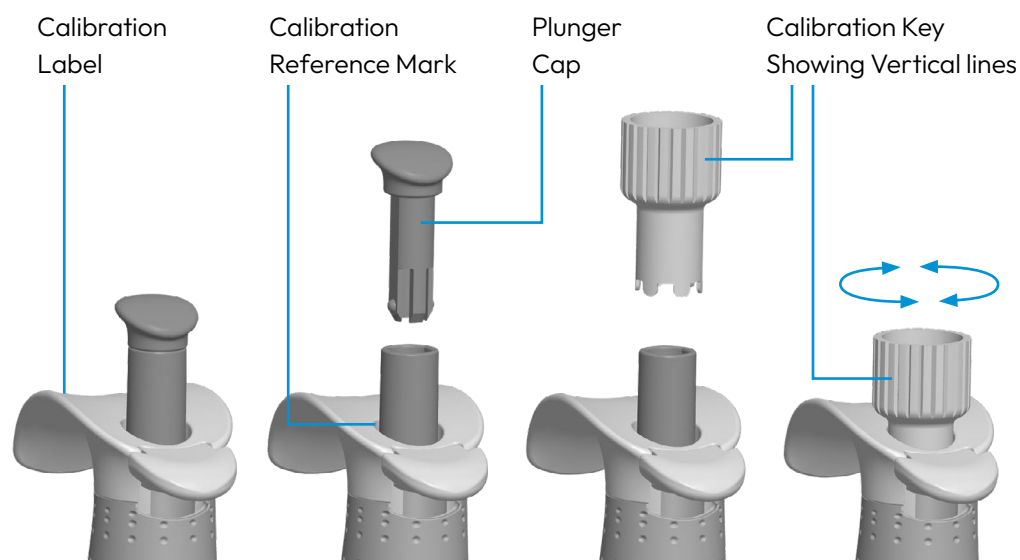

Test Conditions

Liquid	Deionized, degassed, distilled water
Reference Temperature	20-25°C, constant to $\pm 0.5^\circ\text{C}$
Number of Measurements	10 per channel per volume setting, referencing ASTM 1154 Guidelines
Pipette Tips	Hamilton pipette tips

Performance Specifications

Adjustable Volume Pipettes			At 10% of Pipette Volume		At 100% of Pipette Volume	
Part No.	Volume Range	Color Coding	Accuracy within $\pm \%$	Precision within $\pm \%$	Accuracy within $\pm \%$	Precision within $\pm \%$
55019-50	5-50 $\mu\text{L} \times 8$	Sandstone	2.00	2.00	0.80	0.35
55019-52	30-300 $\mu\text{L} \times 8$	Brick Red	1.20	0.70	0.50	0.30
55019-54	5-50 $\mu\text{L} \times 12$	Sandstone	5.00	2.00	1.00	0.50
55019-56	30-300 $\mu\text{L} \times 12$	Brick Red	2.40	0.70	1.00	0.50

Optimizing Performance

Topic	Action
Consistent Technique	Pipette with a consistent rhythm, pressure, and speed.
Tip Selection	Use Hamilton pipette tips for guaranteed accuracy. For information, visit www.hamiltoncompany.com/pipettes .
Tip Size and Fit	Use proper size tip, firmly placed on pipette shaft.
Pre-Rinse Tips	Pre-rinse pipette tips for improved precision.
Sample Aspiration	Keep the disposable tip immersed in fluid during aspiration. Do not let the plunger snap back to starting position.
Immersion Depth	Maintain an immersion depth of 2 to 4 mm.
Immersion Angle	Maintain an angle of <20° from vertical.
Viscous Samples	Aspirate slowly. If bubbles are observed, resample. Volume errors may still occur. Refer to the Calibration section.
Acid Samples	Pipetting strong acids and corrosive solutions is not recommended. These liquids may damage the piston and seal. To protect the pipette, use filter barrier pipette tips and inspect piston frequently. For more information, visit www.hamiltoncompany.com/pipettes .
High Vapor Pressure Samples	Pipetting solutions with high vapor pressure is not recommended. These liquids may damage the piston and seal.
Sample Temperature	Fluids at a temperature other than that for which the pipette and pipette tips have been calibrated may result in volume measurement errors.
Storage	Store upright in stand. Do not lay the pipette on its side with fluid in the tip. Fluid reaching the piston causes contamination and possible corrosion.
Cleaning Piston	Wipe piston with alcohol and a soft, lint-free cloth. Dry and lightly lubricate the piston (Lubricant, Part No. 5763-01).
Seal Replacement	Replace the PTFE seals every 3-6 months depending on use and the samples aspirated.
Performance Checks	Check the accuracy and precision of your pipette every 3-6 months depending on use and the samples aspirated.

Troubleshooting

Problem	Possible Cause	Recommended Remedy
Inaccurate Results	Pipette tip not attached correctly.	Check for tip fit and check for proper tip size, replacing tip as needed.
	Aspiration stroke too rapid (viscous fluids).	Aspirate slowly by controlling rate of plunger release.
	Calibration required.	Calibrate the pipette.
	Seal leak (dripping fluid) .	Replace the seal.
Imprecise Results	Inconsistent sampling method.	See section on Optimizing Performance.
	Seal leak.	Replace the seal.


Cleaning

Wipe the outer surface of the pipette with a mild soap and water daily. In case of accidental contamination of the piston chamber shaft, immediately disassemble the piston chamber. Refer to the Replacement Parts section for instructions. In addition, rinse the ejector plate and piston assembly with mild soap and water followed by distilled water. Dry all components completely. Lightly lubricate the pistons (Lubricant, Part No. 5763-01).

Autoclaving

The entire pipette is autoclavable at 121°C for 15 minutes at 15 psi (103.5 kPa or 1.02 atm) pressure without the need for any disassembly. Although autoclaving should not affect pipette accuracy, practice good laboratory procedures by checking your pipette and recalibrating as needed.

Pipette Parts Diagrams

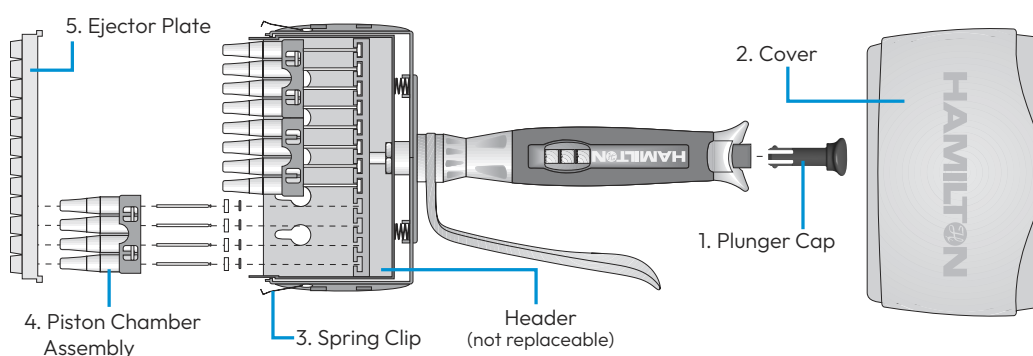

Replacement Parts Table

	Part Description	8 x 50 μ L	12 x 50 μ L	8 x 300 μ L	12 x 300 μ L
1.	Plunger Cap	54190-06	54190-06	54190-10	54190-10
2.	Cover	4917-01	4916-01	4917-01	4916-01
3.	Spring Clip	34354	34354	34354	34354
4.	Piston Chamber Assembly	6385-01	6385-01	6385-02	6385-02
5.	Ejector Plate	4923-01	4922-01	4923-01	4922-01

Replacement Parts

Hamilton pipettes have been uniquely designed so simple repairs and part replacements can be performed by the user. The Replacement Parts Table lists all pipette replacement parts identified by Hamilton part numbers. To replace parts refer to the Pipette Parts Diagram.

1. Snap the front and back covers off of the pipette. If necessary, insert a flathead screwdriver into slot on the side of the lower body and rotate gently.
2. Unhook the two spring clips securing the ejector plate.
3. Remove the ejector plate.
4. Using a Phillips head screwdriver, remove the screw that secures the piston chamber assembly and gently remove piston chamber assembly.
5. Lightly lubricate entire circumference of pistons using a soft, lint-free cloth (Lubricant Part No. 5763-01).
6. Insert and slide the pistons into the new piston chamber assembly. Make sure that the seal flanges are oriented towards the plunger with the O-rings below the flanges towards the tip.
7. Adjust the position of the piston chamber assembly so that the screw on the back goes into the large end of the opening on the opposite side. With the screw side facing you, slide the piston chamber down so the screw goes to the very bottom of the opening, and tighten the screw.
8. Reinstall the ejector plate, with the longer lip facing the front (the side without the screws), and secure the ejector plate with the spring clips.
9. Snap the front and back covers back together.

To assure that your instrument meets the original Hamilton Company standards and specifications, we strongly suggest that you follow Good Laboratory Practices by recalibrating your pipette after replacing any parts.

Pipette Recertification and Repairs

Recertification

Hamilton Company offers a program to recertify the accuracy and precision of your Hamilton pipette. Testing is performed at a Hamilton facility under stringent conditions on state-of-the-art equipment. The calibration process is performed with an unbroken chain of calibrations traceable to N.I.S.T. Your pipette will be returned to you with a new Certificate of Performance. If a pipette must be repaired to bring it into conformance, you will be contacted with a cost estimate.

Repairs

Hamilton Company offers a pipette repair program in addition to its general, 12-month warranty. You may return your Hamilton pipette for evaluation. A series of basic tests will be performed to evaluate the pipette's performance. If repairs are required to assure and confirm that your pipette is functioning to original Hamilton specifications, you will be contacted with a cost estimate and the additional repairs and recalibration will be performed upon your approval. Your pipette will be returned to you with a new Certificate of Performance.

For pipette recertification and repairs in the U.S.A., call 1-888-525-2123. For authorized service outside of the U.S.A., contact your local Hamilton Company office or representative using the contact information on the back cover.

Return of Goods

Hamilton return and repair policy is written to protect our employees from potentially hazardous materials (e.g., infectious agents, radioactive materials, and carcinogenic chemicals) or any substance that may cause them partial or permanent disability during the inspection or repair process. When returning products, please ensure the items are decontaminated and marked accordingly. Furthermore, the customer assumes responsibility should the returned product be determined to be hazardous. All returns must be accompanied by a pre-approved Returned Goods Authorization Number (RGA#). Any return not received without a RGA# may be disposed. Contact Hamilton Customer Service at 1-888-525-2123 for an RGA#.

Hamilton Precision Pipette Tips

Hamilton delivers performance, value, and choice with two high quality tip families, plus multiple formats and packaging options. The tip families are low retention/low binding tips for maximum sample recovery and standard tips for routine, high-throughput work. Formats include standard, non-filter, extended length, hydrophobic filter barrier, non-sterile, pre-sterilized, and macrovolume. Packaging options include bulk bags, hinged and lift-off racks, and stackable racks. Hamilton tips are made from chemically resistant, virgin polypropylene for excellent clarity and smooth surfaces. All Hamilton Pipette tips are certified to be free of RNase, DNase, and pyrogens. Hamilton pipette tips are universal-fit. For more information, visit www.hamiltoncompany.com/pipettes.

The following are trademarks of Hamilton Company:

AdvanTip™
Pipette Safe™
SoftGrip™
SofTouch™
SofTop™
SoftAide™

Quality Hamilton Products:

MICROLITER™ Syringes
GASTIGHT® Syringes
Chromatography Syringes
Syringes for Life Science
Instrument Syringes
SoftGrip™ Pipettes
Miniature Valves
Modular Valve Positioner
Laboratory Fittings,
Adapters & Tubing
Precision Syringe Pumps
MICROLAB® Diluters & Dispensers
Electrochemical Sensors
DURACAL™ Buffer Solutions
HPLC Columns & Resins

Laboratory Automation for:

Drug Discovery
Genomics
Proteomics
Forensics
In Vitro Diagnostics



● Headquarters / Manufacturing



Years of Experience
75+



Locations Worldwide
22+



Employees Internationally
3,000+

To find a representative in
your area, please visit:
www.hamiltoncompany.com/contacts

Web: www.hamiltoncompany.com
USA: 800-648-5950
Europe: +40-356-635-055

Hamilton Americas & Pacific Rim
Hamilton Company Inc.
4970 Energy Way
Reno, Nevada 89502 USA
Tel: +1-775-858-3000
Fax: +1-775-856-7259
sales@hamiltoncompany.com

Hamilton Europe, Asia & Africa
Hamilton Central Europe S.R.L.
str. Hamilton no. 2-4
307210 Giarmata, Romania
Tel: +40-356-635-055
contact.hce.ro@hamilton-ce.com

